



REVISIONS			
ZONE	REV	DESCRIPTION	DATE
		APPROVED	

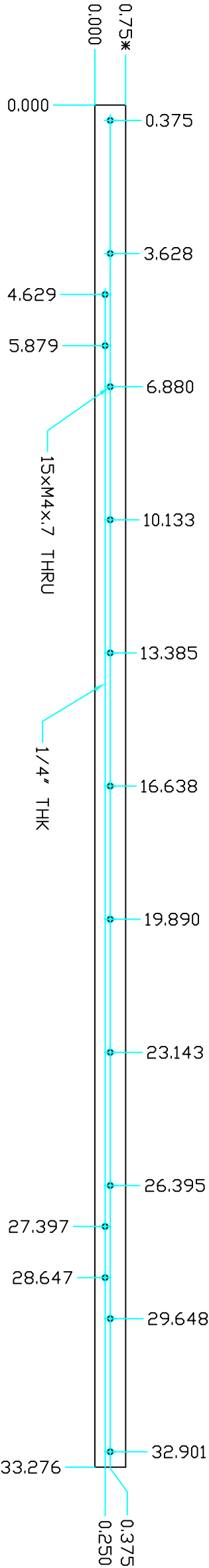

PART 1

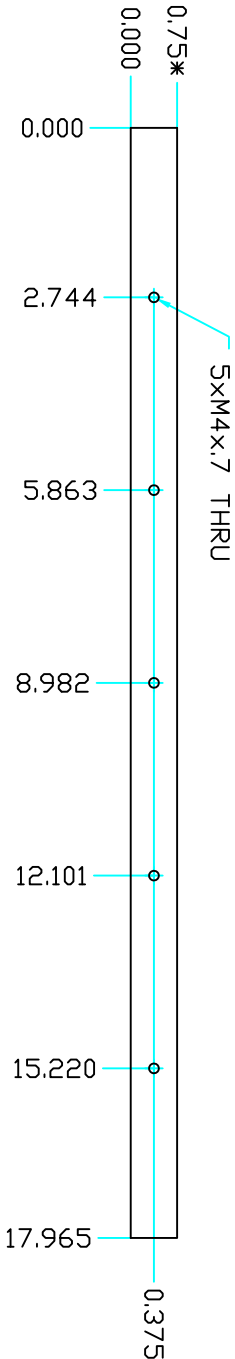

PART 2

2	2	EA	PART 2	ALUMINUM 6061	ALUMINUM BAR 1/4" THK x 3/4" W	103.00.02.02.02
1	2	EA	PART 1	ALUMINUM 6061	ALUMINUM BAR 1/4" THK x 3/4" W	103.00.02.02.01
FIND/REF #		QTY.	UOM	PART NAME	MATERIAL	REF. SHEET/DWG.
PARTS LIST						
UNLESS OTHERWISE SPECIFIED				PERMISSIBLE VARIATION IN INCH DIMENSIONS WITHOUT TOLERANCES		
1. DIMENSIONS ARE IN INCHES.				FABRICATION TOLERANCES		
2. MACHINE SURFACES TO BE FLAT, SQUARE, PARALLEL, & CONCENTRIC BELOW 12				±.010	BETWEEN PLACES X & X.X	X.XXX
3. WITHIN .001 FOR .5 INCHES OR .001 TIR AS APPLICABLE.				12 TO 60	±.013	±.010
4. HOLE LOCATIONS (EXCEPT COMMERCIAL PARTS) WITHIN .001				60 TO 180	±.020	±.010
5. BREAK ALL SHARP CORNERS .01 x 45° CHAMFER.				OVER 180	±.030	±.020
6. FILLETS SHOWN SHARP = .04 RAD.				OVER 48	±.045	±.010
7. TAPPED HOLES UNITED CLASS 2B THREAD, MIN DEPTH 1.5 x DIA.				ANGLE TOLERANCE	±.01°	
8. CENTER PERMISSIBLE						
9. HOLE LOCATIONS WITHIN .001						
10. ALL WELDING TO CONFORM TO LATEST AWS STANDARDS						
SURFACE TEXTURE				125		
TOLERANCES ARE NOT ACCUMULATIVE						

103.00.02.02.02

SIZE	SCALE
D	1:4
PROJECT NUMBER	
103	
DRAWING NUMBER	
103.00.02.02	
REVISION	
1 OF 1	

Open Source Instruments Inc.

FARADAY ENCLOSURE, FE33

OSI

\* REFERENCE ONLY

