


REVISIONS			
ZONE	REV	DESCRIPTION	DATE
APPROVED			

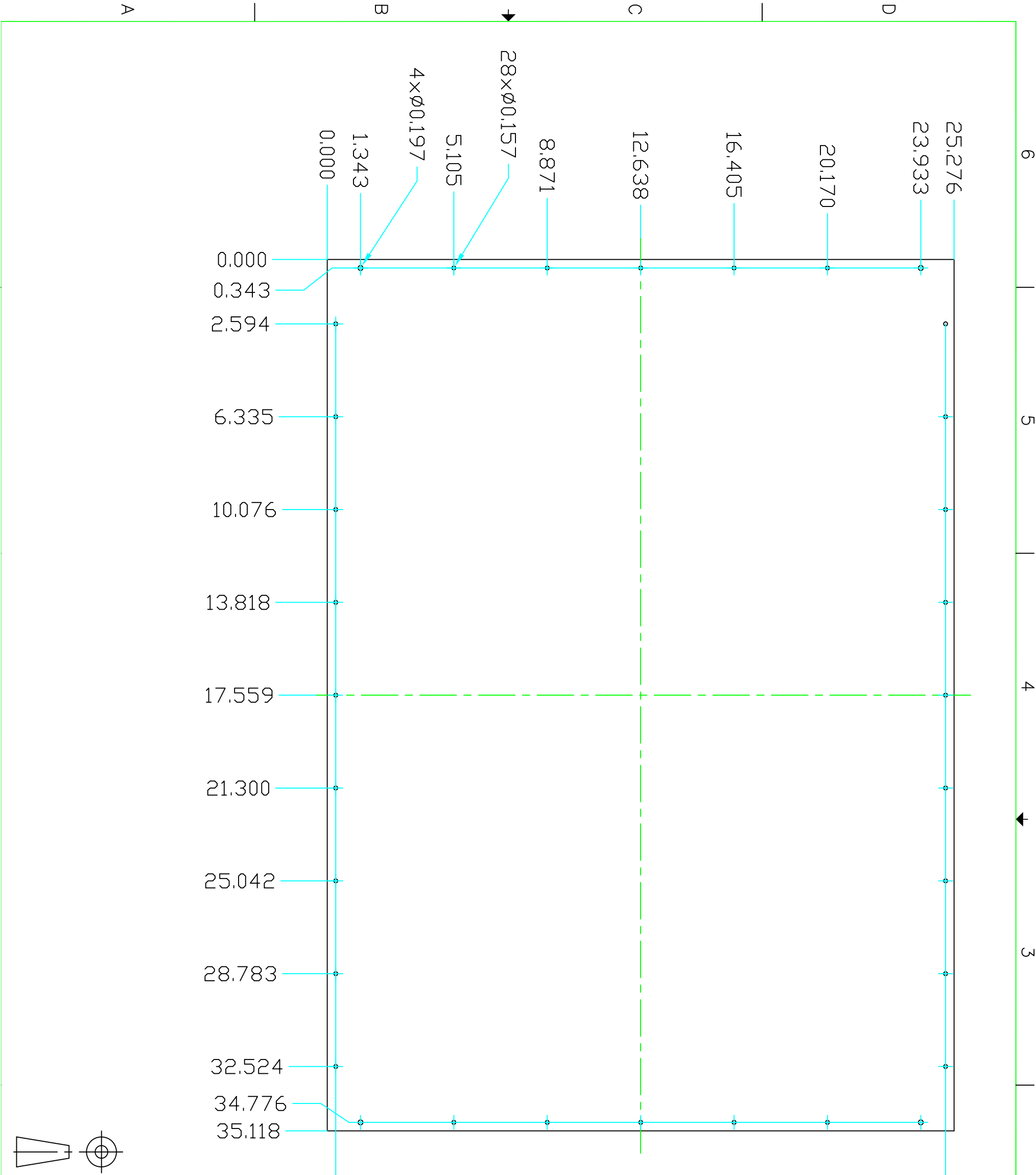

1	2	EA	PART 1	ALUMINIUM 6061 24"x24"x0.032 THK SHEET	102.00.01.01
FINI/REF.#	QTY.	UM.	PART NAME	MATERIAL	REF. SHEET/DWG.
PARTS LIST					
UNLESS OTHERWISE SPECIFIED				PERMISSIBLE VARIATION IN INCH DIMENSIONS WITHOUT TOLERANCES	
1. DIMENSIONS ARE IN INCHES				FABRICATION TOLERANCES	
2. MACHINE SURFACES TO BE FLAT, SQUARE, PARALLEL, & CONCENTRIC				MACHINING TOLERANCES	
3. MARK ALL DETAILS (EXCEPT COMMERCIAL PARTS) WITH DWG.				BELOW 12	
4. BREAK ALL SHARP CORNERS .01" X 45° CHAMFER				12 TO 60	
5. TOLERANCES SHOWN SHARP = .04 RAD				60 TO 180	
6. TOLERANCES SHOWN SHARP = .04 RAD				18 TO 48	
7. CENTER HOLE DIA. 1/16" X .03"				OVER 48	
8. CENTERS PERMISSIBLE				ANGLE TOLERANCE ±0.1°	
9. DO NOT SCALE DRAWING				SURFACE TEXTURE: 125	
10. ALL WELDING TO CONFORM TO LATEST AWS STANDARDS				TOLERANCES ARE NOT ACCUMULATIVE	
DRAWING INFORMATION				SIZE: D	
NAME: AD				SCALE: 1:5	
CHECKED: KH				PROJECT NUMBER: 103	
APPROVED: KH				DRAWING NUMBER: 103.00.00.01	
TITLE: FARADAY ENCLOSURE, FEAS3 WALL					
OSI					
SHEET: 1 OF 1					
REVISION: A					