
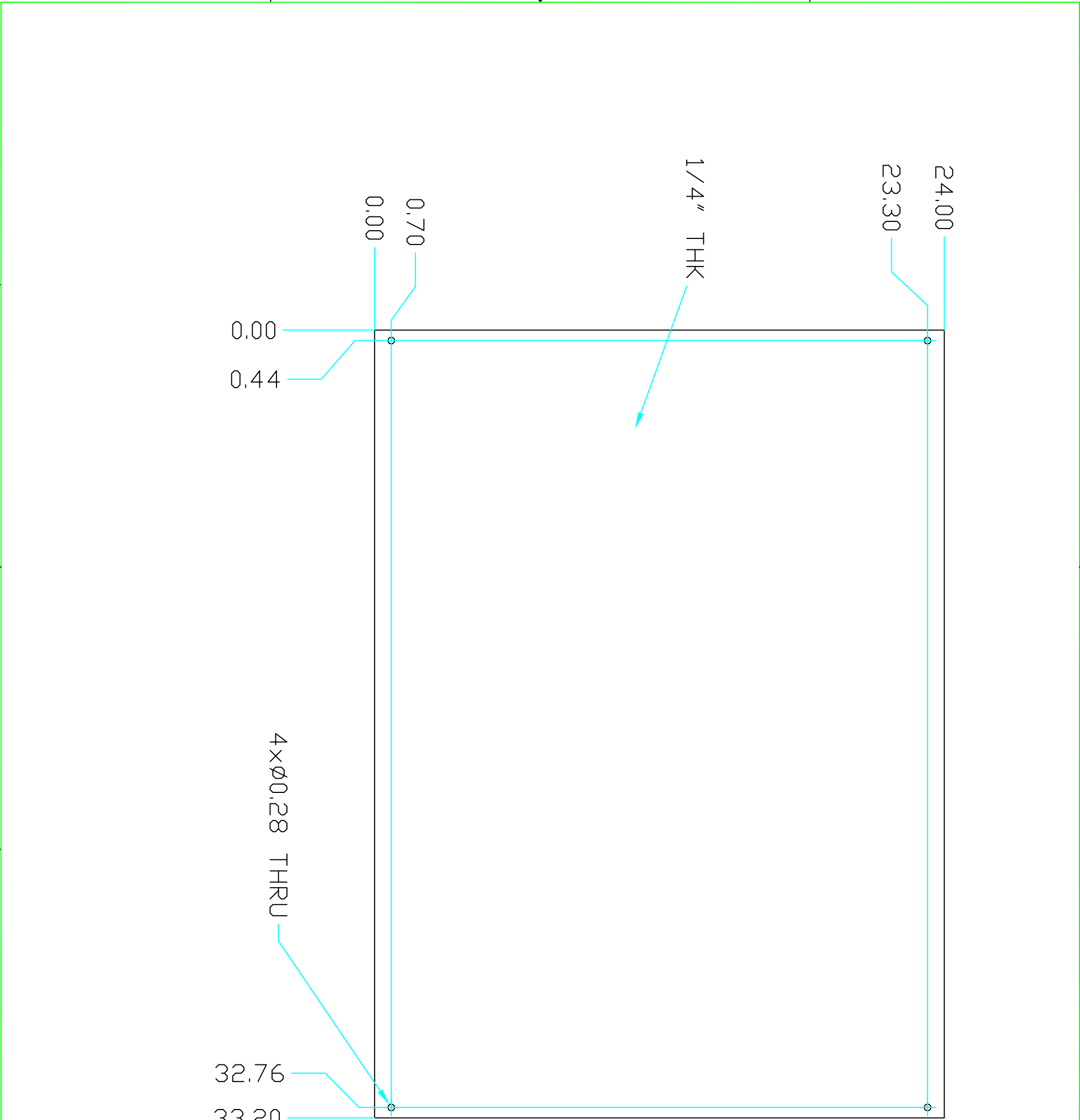


REVISIONS		DATE	APPROVED
ZONE	REV	DESCRIPTION	

END/REF#	QTY.	U.M.	PART NAME	MATERIAL	CLEAR POLYCARBONATE SHEET 36" L x 24" W x .023" THK	REF. SHEET/DWG.			
PARTS LIST									
UNLESS OTHERWISE SPECIFIED									
FABRICATION TOLERANCES									
MACHINING TOLERANCES									
PERMISSIBLE VARIATION IN INCH DIMENSIONS WITHOUT TOLERANCES									
1. DIMENSIONS ARE IN INCHES									
2. MACHINING SURFACES TO BE FLAT, SQUARE, PARALLEL & CONCENTRIC, BELLY 12									
3. MARK ALL DETAILS (EXCEPT COMMERCIAL PARTS) WITH DWG.									
4. BREAK ALL SHARP CORNERS .01 x 45° CHAMFER									
5. FILET'S SHOWN SHARP = .124 RAD									
6. FINISH ALL SURFACES TO CLASS 25 THREAD, MIN DEPTH .15 x DIA									
7. CENTER HOLE & REAR HOLE 1/16 x 45°									
8. CENTERS PERMISSIBLE									
9. DO NOT SCALE DRAWING									
10. ALL WELDING TO CONFORM TO LATEST AWS. STANDARDS									
SURFACE TEXTURE 125									
TOLERANCES ARE NOT ACCUMULATIVE									
DRAWING INFORMATION									
NEXT ASSY	QTY	NAME	DATE	Open Source Instruments Inc.			SIZE	SCALE	
		BRN	AD	06/22/23				D	1:4
		CHECKED	KH					PROJECT NUMBER	
		APPROVED	KH					109	
TITLE							DRAWING NUMBER		
FARADAY ENCLOSURE, FE3B							109.00.00.06		
SHEET							REVISION		
1 OF 1							A		

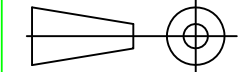

6

5

4

3

2

1